



Prof. Dr. R. Verena Taudte
Telefon: +49 6421 58-66953
Dr. Jelena Pesek
Telefon: +49 6421 58-64096
Lena Ludwig-Radtke
Telefon: +49 6421 58-67018
E-Mail: metabolomics@uni-marburg.de



Vorgaben für die Vorbereitung von Proben zur massenspektrometrischen Analyse in der Core Facility für Metabolomics

- Probengefäße müssen resistent gegenüber -80°C sein
- Bitte nur folgende Probengefäße verwenden:
 - o Reaktionsgefäße („Eppis“) mit Safe-Lock-Deckel
 - o Probenfläschchen (GC-Vials) Volumen 1,5 ml; Durchmesser 12 mm
- Die Beschriftung muss mit Etiketten oder lösungsmittelresistenten Stiften erfolgen
- Proben dürfen ausschließlich fortlaufend nummeriert sein (Deckel und Seite)
- Bitte keinen Parafilm verwenden
- Die Proben müssen in beschrifteten Boxen (nicht in Tüten) einsortiert werden
- Bitte vor jeder Probenabgabe eine aktuelle Probenliste senden

Guidelines for preparing samples for mass spectrometry analysis in the Metabolomics Core Facility

- Sample containers must be suitable for -80°C storage
- Please use only the following sample containers:
 - o Reaction tubes („eppis“) with safe-lock lids
 - o Sample vials (GC vials): volume 1.5 ml; diameter 12 mm
- Labelling must be done with labels or solvent-resistant markers
- Samples may only be numbered (lid and side)
- Do not use Parafilm
- Samples must be sorted in labelled boxes (not in bags)
- Please provide an updated sample list before each sample submission